FACT SHEET

HEALTH, SAFETY, AND RISK MANAGEMENT

UNIVERSITY OF MINNESOTA Driven to Discover®

Proper Removal of Contaminated Gloves



Figure 1: Grip one glove near your wrist.

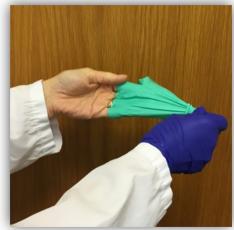


Figure 3: Pull the glove off & grip it in your gloved hand. Do not snap the glove.



Figure 5: Pull the 2nd glove off the tips of your fingers, concealing the 1st glove inside.



Figure 2: As you pull the glove over your fingers, the glove will be pulled inside-out.



Figure 4: Slide your finger under your 2nd glove, still gripping the 1st glove.



Figure 6: The 1st glove will be contained within the 2nd glove. Dispose appropriately.